IKAMAG® RET basic
safety control

IKAMAG® RET basic C
safety control

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OPERATING INSTRUCTIONS EN 9

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DE

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CE-DECLARATION OF CONFORMITY

EN

We declare under our sole responsibility that this product corresponds to the regulations 89/336/EG and 73/23/EG and conforms with the standards or standardized documents DIN EN IEC 61010-1 and DIN EN IEC 61326-1.

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Warranty

You have purchased an laboratory machine which meets the highest engineering and quality standards.

In accordance with guarantee conditions, the guarantee period is 24 months. For claims under the guarantee please contact your local dealer. You may also send the machine direct to our works, enclosing the delivery invoice and giving reasons for the claim. You will be liable for freight costs.

The warranty does not cover wearing parts, nor does it apply to faults resulting from improper use or insufficient care and maintenance contrary to the instructions in this operating manual.

If you require servicing, return the appliance in its original packaging. Storage packaging is not sufficient. Please also use suitable transport packaging.

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Contens

	Page
Warranty	2
Safety instructions	9
Proper use	10
Unpacking	10
Commisioning	10
Switching on	10
Function Heating	10
Adjusting the safety temperature limit "SAFE TEMP"	11
Control of the average temperature with the elec- tronic temperature sensor	11
Function Stirring	12
Maintenance and Cleaning	13
Associated standards and regulations	13
Accessories	13
Technical data	14
Information for care and maintenance of the heating plate vers. C	21
List of spare parts RET basic/C	22
Spare parts diagram RET basic/C	23

Safety instructions

Caution: This device must only be used to process and heat media whose flash point is over the safe temperature limit set point (50 to 380°C).

The adjusted safety temperature must always be appropriate for 25°C below the flash point of the used liquid.

Examine regularly (we recommend every 4 weeks) the function of the safety temperature limiter. For this the safety temperature must be changed around at least 2 scale parts of the value already stopped and be accomplished the attitude as under

“adjusting the safety temperature delimitation SAFE temp” decribed. By the exceeding of the adjusted safety temperature (e.g. rotary button “temp” posed on max.) you can examine the function. The “temp” LED must expire and the heating must be lasting switched off.

The instrument may not be operated in rooms with explosion hazard.

Furthermore, it has to be placed upon a fire-proof and/or not burnable surface.

Attention - Magnetism! Effects of the magnetic field have to be taken into account (e.g. data carriers, cardiac pacemakers ...).

A separation from the line is made with the equipment only by pulling net and /or device plugs. The plug socket for mains connection line must be easily attainable and accessible.

Please make sure that the mains cable does not contact the heating plate. The power cable must be replaced by a cable of equal quality and capacity.

When using PTFE-coated magnetic bars, the following hase to be noted: *Chemical reactions of PTFE occur in contact with molten or dissolved alkaline and alkaline - earth metals, as well as with fine-particled powders of metals of the 2. and 3. group of the peridocal system at temperatures above 300-400°C.*

Only elementary fluorine, chlorine trifluoride und alkaline metals do attack PTFE, halogen hydrocarbons have a reversibly swelling effect.

Source: Römpps Chemie-Lexikon and „Ullmann“ Bd.19

Exercise caution when touching the housing parts and the heating plate!

Risk of burns!

The heating plate can reach temperatures in excess of 340°C.

Please consider possible endangering due to release of toxid or flammable gaser caused by heated medium.

Please note that wear debris from the appliance or rotating accessories can end up in the medium to be processed. Please contact **IKA** should you have any questions in this respect.

Proper use

The RET basic or RET basic C is a stirring instrument with heating function. Combined with the ETS-D electronic temperature sensor, the RET basic or RET basic C can be used as a **safety magnetic stirrer**.

It is used in e.g. laboratories of the chemical industry, in schools as well as in pharmacies. The instrument is suitable for tempering substances which are placed on the heating plate in vessels. By means of a built-in stirring drive and with the help of a magnetic bar within the vessel, the substances can be stirred at the same time. The mixing intensity is dependent on the speed and on the size of the magnetic bar.

Unpacking

Please unpack the equipment carefully and check for any damages. It is important that any damages which may have arisen during transport are ascertained when unpacking. If applicable a fact report must be set immediately (post, rail or forwarder).

The delivery scope covers: an RET basic or RET basic C, connecting cable and operating instructions.

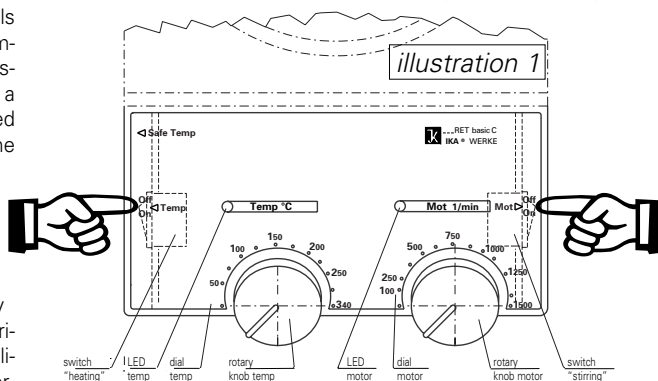
Commissioning

Check whether the voltage specified on the type plate matches the mains voltage available. The power socket used must be earthed (protective earth conductor contact). If these conditions are met, the device is ready to operate after plugging in the mains plug. If these procedures are not followed, safe operation cannot be guaranteed and/or the equipment may be damaged.

Observe the ambient conditions (temperature, humidity, etc.) listed under Technical Data.

Switching on

After the instrument has been switched on, it is in „standby“-operation. The functions „heating“ and „stirring“ are started with the help of switches on the left and right bottom of the housing (see illustration 1)

Function Heating

The heating plate temperature of the instrument is kept at a constant temperature by the control circuit. The heating plate temperature is also monitored by a separate adjustable safety circuit. The temperature sensors required for this (two Pt 100's) are integrated into the heating plate.

The heating plate has an output of 600 Watts.

The LED next to the imprint "Temp" will shine green when turning on the function "heating". The temperature of the heating plate can be adjusted at the rotary knob "Temp" with the included scale (0 up to 340°C). During the heating period (phase) the colour of the LED "Temp" will change from green into orange (orange = supp-

ly of energy to the heating plate, green = the adjusted temperature is achieved). Dependent on the required heating capacity, the "Temp"-LED can blink in a 0,5 second interval.

If there is a malfunction in the control circuit, the heating plate will be **permanently tuned off** by the safety circuit.

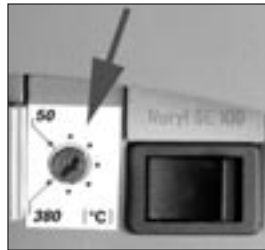
PLEASE NOTE:

The heating plate will not heat up unless the contact plug is plugged in (the contact plug is permanently attached to the instrument) or a **ETS-D** or **contact thermometer** is connected.

Adjusting the safety temperature limit "Safe Temp"

ACHTUNG: The adjusted temperature must always be appropriate for 25°C below the flash point of the used liquid.

The "Safe Temp" set screw is located on the left next to the "Temp" adjustment dial under the control panel.

The safety temperature (max. heating plate temperature) can be adjusted with the "Safe Temp" set screw of the safety temperature limiter. As soon as the desired safety temperature is exceeded by a higher adjusted set-point temperature or a defect

occurs in the temperature controller, the "Temp" LED goes out. The heater turns off permanently in this status and **the stirring function continues to run at the speed that was set before the malfunction.**

This adjustment has to be carried out without a vessel on the heating plate.

- Switching on the heating with the rocker switch "OnOff/Temp"
 - Turn the "Safe Temp" set screw as far to the right (Clockwise) as possible with a straight-blade screwdriver (=ca. 380°C).
 - Adjust the set point temperature to the desired safety temperature (Safe "Temp" with the "Temp" dial).
 - Adjust the set point temperature to the desired safety temperature ("Safe Temp") with the "Temp" dial.
Example: Of a liquid with a flash point when using 125°C the target temperature (rotary button TEMP) may be stopped maximally to 100°C.
 - Please wait till the heating reaches the requested temperature (safety temperature) (this may take some minutes), the "Temp"-LED will blink (approx. 2-times per second).
 - Turn the "Safe Temp" set screw slowly to the left up to the point where it switches off. Then and only then, the green "Temp" LED goes out.
 - Let the heating plate cool off.
 - After you have turned the instrument off and back on again with the "Temp On-Off" switch, it is ready for operation.
- Note:** If the "Safe Temp" set screw is set to minimum (as far as will go to the left), the "Temp" LED goes out and the heater is turned off).

Control of the average temperature with the electronic temperature sensor

The preferable method for controlling the average temperature is with the ETS-D or contact thermometer. After the set point temperature has been adjusted, this results in a short heating-up time, practically no temperature drift and only minor fluctuation in temperature.

In addition to precise control function, the adjustabel "MAX-TEMP" function makes the ETS-D a temperature limiter that has a direct effect on the average temperature. When this "MAXTEMP" is exceeded, for example because of a defect in the control circuit or unintentional of the "Safe Temp" or set

point temperature on the RET basic, the ETS-D immediately and permanently turns off the separate circuit of the RET basic.

The stirring function then continues to run at the speed that was set before the malfunction.

A 6-pin jack is located on the rear side of the instrument for connecting the ETS-D, the contact thermometer or the contact plug. The electronics of the instruments returns a test current that must flow via connector pins 3 and 5 for the heating plate to heat up.

Safety contact thermometers

acc. to DIN 12 878 class 2 or acc. Gerstel are connected with a three-wire cable, the test current flows through the contact thermometer.

Safety function:

If the test current is interrupted because of e.g. breakage of contact thermometer or falling out of the cable plug, the heating cuts off. The required plug allocation is indicated on a reference plate above the diode bushing.

Contact thermometer without safety circuit

acc. to DIN 12 878 class 0. The instrument only heats if the test current circuit is closed by an electrical connection of the plug pins 3 and 5.

2-wire connecting cables:

Connect plug pins 3 and 5 of the instrument plug.

3-wire connecting cables:

Here the test current circuit can also be produced in the terminal head of the contact thermometer (connect plug pins 2 and 3). A 3-wire cable with the required bridge is available (accessories).

Settings

For the detailed instructions for settings and limit values, please refer to the operating instructions for the instrument you are connecting.

The desired average temperature can be adjusted on the ETS-D or contact thermometer. In addition, a "MAXTEMP" can be set for the ETS-D, in other words, a maximum average temperature can be set for malfunctions in combination with the RET basic. The required surface temperature of the heating plate can be selected with the "Temp" dial.

Adjusting the "Temp" dial to the maximum adjustable temperature will result in the fastest possible heating time, but the average temperature may fluctuate to values above the set-point temperature on the contact thermometer, for example. By adjusting the "Temp" dial to approximately twice the set-point value (thus with a set-point of 60°C the Temp dial would be set to 120°C), you will reach a good compromise between a fast heating time and overshooting the set point.

If you adjust the Temp dial to exactly the set-point temperature, the medium will not reach the set-point temperature, since some loss of the heat will always occur between the heating plate and the medium.

The maximum heating plate temperature can be adjusted with the "Safe Temp" screw.

Function Stirring

The function stirring is started and stopped by its corresponding switch. If the function "Stirring" is started, the LED "Mot" beside the the imprint "Mot" shines green.

The speed of the drive can be adjusted on the "Mot" dial with the corresponding scale from 50 to 1500 rpms.

The motor speed of the instrument is kept constant by a control circuit.

Maintenance and Cleaning

The RET basic and RET basic C is maintenance-free. It is subject only to the natural wear and tear of components and their statistical failure rate.

When ordering spare parts, please give the manufacturing number shown on the type plate, the machine type and the name of the spare part.

Please send in equipment for repair only after it has been cleaned and is free from any materials which may constitute a health hazard. Use only cleansing agents which have been approved by IKA to clean IKA devices. To remove use:

Dyes	isopropyl alcohol
Construction materials	water containing tenside / isopropyl alcohol
Cosmetics	water containing tenside / isopropyl alcohol
Foodstuffs	water containing tenside
Fuels	water containing tenside

For materials which are not listed, please request information from IKA. Wear the proper protective gloves during cleaning of the devices.

Electrical devices may not be placed in the cleansing agent for the purpose of cleaning.

Before using another than the recommended method for cleaning or decontamination, the user must ascertain with the manufacturer that this method does not destroy the instrument.

Protective conductor:

On the under side of the heating plate are areas (bare metal) and can be used as a test point for the protective conductor. It may be necessary to remove an oxide coating for the subsequent tests.

Associated standards and regulations

Construction in accordance with the following safety standards

EN 61010-1	EN 61010-2-10
UL 3101-1	CAN/CSA C22.2 (1010-1)

Construction in accordance with the following EMC standards

EN 61326-1+A1

Associated EU guidelines

EMC-guidelines: 89/336/EG

Machine guidelines: 73/23/EG

Accessories

RS 1	set of stirring bars	H16V	support rod
RSE	PTFE stirring bar remover	H 15	bath top
H 44	cross sleeve	TC 1	Electronic contactthermoter
H 36	holding rod		

Recommended IKA-stirrers

TRIKA-stirrers PTFE, triangular
lengths 25, 40 and 55 mm

IKAFLON-stirrers PTFE
dia 6mm lengths 10 and 15 mm
dia 8mm lengths 20, 25, 30, 40, 50 and 60 mm

Technical data

design voltage:	VAC 220 - 230 ±10%
oder	VAC 115 ±10%
oder	VAC 100 ±10%
design frequency:	Hz 50/60
input power total	W 630 (by 230V)
Drive:	EC-motor with ball bearings, without brushes
Max. power consumption / output:	W 12 / 5
speed range:	rpm 50...1500 infinitely variable setting
speed variety (without load, 1500rpm):	% ± 10
stirring quantity (H ₂ O):	l 20
Heating plate:	
diameter of heating plate:	mm 135
heating power: (230V / 115V)	W 600 ±10%
(100V)	W 455 ±10%
temperature range:	°C RT 340 ±20K
Infinitely variable safety temperature "SAFE TEMP"	°C 50 - 380 ±20K
temperature fluctuation of the surface temperature	K ±1
control accuracy of medium in ETS-D operation	K ±1
contact thermometer- connection	DIN 12 878 class 2
permitted ambient temperature:	°C +5 to +40
permitted relative humidity:	% 80
permitted duration of operation:	% 100
prozection class acc. EN 60529:	IP 42

fuses:	F1/F2 T6,3A /250V Id-Nr 14 080 00 F3 T0,8T /250V Id-Nr 27 562 00
contamination level:	2
overvoltage category:	II
protection class:	1(protective earth)
operation at a terrestrial altitude	m max 2000 above sea level
dimensions:	mm 160 x 280 x 85
weight:	kg 2,4

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Information for Care and Maintenance of the Heating Plate with Technical Enamel Coating

The technical enamel coating makes the heating plate easier to care for and more resistant to acids and bases. Because of it, however, the heating plate is also more susceptible to extreme fluctuations in temperature and the force of impact. This can result in cracks forming or the coating flaking off.

Make certain that the bottom of the placing vessel is even, clean and dry. The bottom of the placing vessel must not have any sharp grooves, sides or edges. Remove residues of bases and acids immediately. We recommend most strongly that you clean the heating plate regularly.

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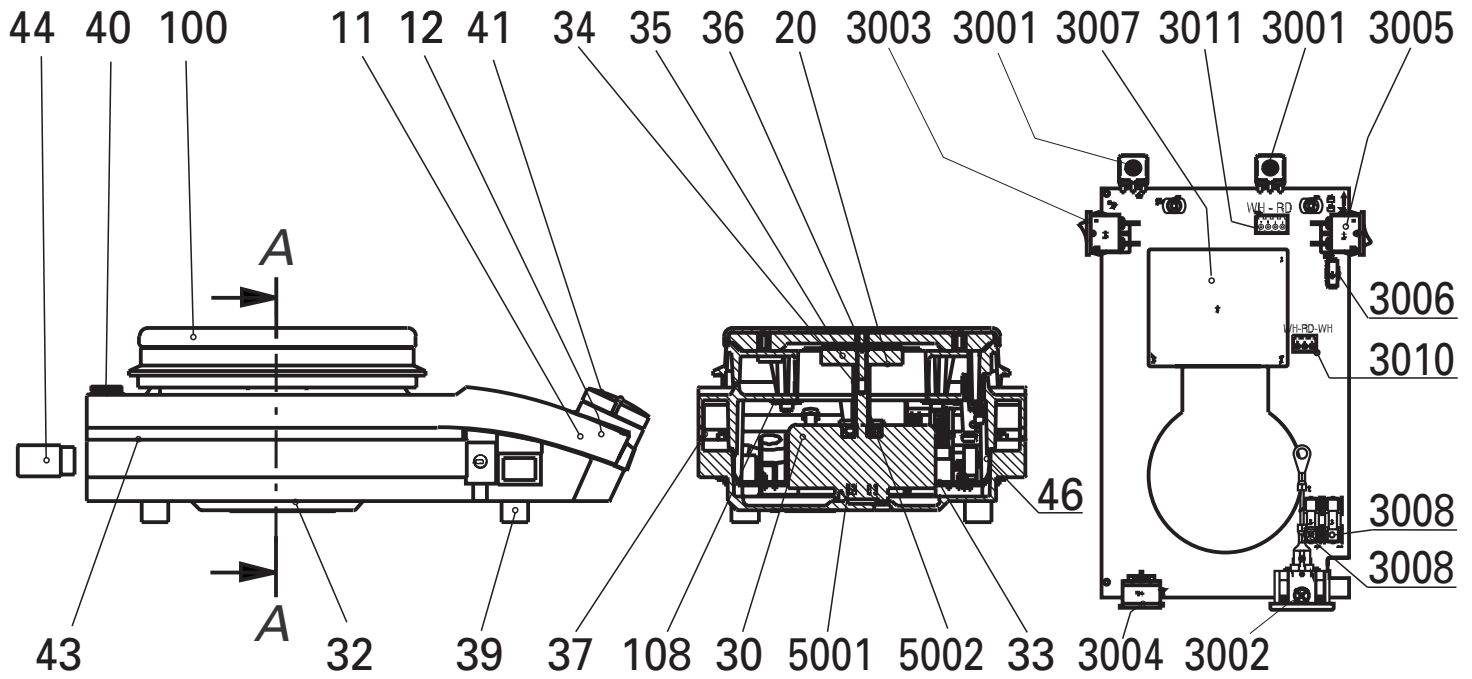
IKAWorks, Inc.

LABORATORY TECHNOLOGY
ANALYZING TECHNOLOGY
PROCESSING EQUIPMENT

North America

IKAWorks, Inc.

2635 North Chase Pkwy SE
Wilmington, NC 28405-7419
USA
TEL. +1 800 733-3037
TEL. +1 910 452-7059
FAX +1 910 452-7693
E-mail: usa@ika.net

IKAWorks, (Asia) Sdn Bhd

LABORATORY TECHNOLOGY
ANALYZING TECHNOLOGY
PROCESSING EQUIPMENT

Asia - Australia

IKAWorks (Asia) Sdn Bhd

No. 17 & 19, Jalan PJU 3/50
Sunway Damansara Technology Park
47810 Petaling Jaya
Selangor, Malaysia
TEL. +60 3 7804-3322
FAX +60 3 7804-8940
E-mail: sales@ika.com.my

IKAJapan Y.K.

LABORATORY TECHNOLOGY
ANALYZING TECHNOLOGY
PROCESSING EQUIPMENT

Japan

IKAJapan Y.K.

293-1 Kobayashi-cho
Yamato Koriyama Shi
639-1026 Japan
TEL. +81 74358-4611
FAX +81 74358-4612
E-mail: japan@ika.de

IKAWorks do Brasil Ltda.

LABORATORY TECHNOLOGY
ANALYZING TECHNOLOGY
PROCESSING EQUIPMENT

South America

IKAWorks do Brasil Ltda.

Estrada do Guerenguê, 491
Taquara Jacarepagua, RJ
Rio de Janeiro
22713-000 Brasil
TEL. +55 21 2435-9600
FAX +55 21 2435-9601
E-mail: brasil@ika.net

IKAWorks Guangzhou

LABORATORY TECHNOLOGY
ANALYZING TECHNOLOGY
PROCESSING EQUIPMENT

China

IKAWorks Guangzhou

173-175 Friendship Road
Guangzhou Economic & Technological
Development Zone
Guangzhou 510730, P.R.CHINA
TEL. +86 20 8222-6772
FAX +86 20 8222-6776
E-mail: sales@ikagz.com.cn