


## IKA® HG 600 IKA® HG 1000

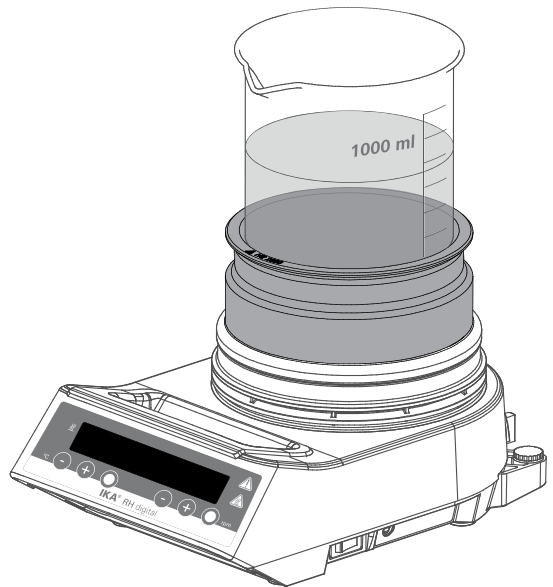

Operating instructions

## Safety instructions



• Read the operating instructions in full before starting up and follow the safety instructions.

• Keep the operating instructions in a place where it can be accessed by everyone.  
• Follow the safety instructions, guidelines, occupational health and safety and accident prevention regulations.



• Pay attention to the residual heat after processing. The heating jacket temperature may be very high.

• Allow it to cool down when handling the heating jacket.

• Never carry the heating jacket around as long as they are hot.

• The heating jacket and the beaker must be properly placed on the heating plate and cannot come off by themselves. The centre of gravity of the heating jacket and the beaker must lie in the heating plate.

## General information

**IKA® HG 600 or HG 1000** heating jacket use for matching with **IKA®** magnetic stirrer with heating or heating plate. The heating jacket improves heat transfer into the medium. It can cut down the heating time considerable.

Anodizing color of the heating jacket will change with aging, it do not affect normally heating function of the heating jacket.

## Cleaning

Use only cleaning agents which have been approved by **IKA®** to clean the heating jacket.

### Dirt

Dye

Construction material

Cosmetics

Foodstuffs

Fuel

### Cleaning agent

isopropyl alcohol

water containing tenside / isopropyl alcohol

water containing tenside / isopropyl alcohol

water containing tenside

water containing tenside

## Technical data

		HG 600	HG 1000
For beaker volume	ml	600	1000
Material		Aluminium	
Inner diameter	mm	91	106
Dimensions	mm	Ø 111 x 50	Ø 126 x 50
Weight	kg	0.35	0.40

### IKA® - Werke GmbH & Co.KG

Janke & Kunkel-Str. 10


D-79219 Staufen

Tel. +49 7633 831-0

Fax +49 7633 831-98

sales@ika.de

[www.ika.com](http://www.ika.com)

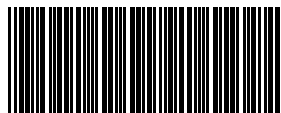

\*25001641\*